

Download and Installation procedure for 30 days free trial version

- ① Download , in any folder, 3 files from the site below;
http://www.yokogawa.com/scanner/product/ca/ca_6_e.htm

Download

[Top](#)
[Workflow](#)
[Application](#)
[Download](#)

■ Download

Software, User's Manual, Tutorial, Protocol guides and Sample data are available for downloading.

Activation code is required to start the software. Please download the requesting form and send it to the e-mail address described on the form.

| Products | Download | | | | | | |
|---------------------------------------|--|-----------------------|-------------------------------|----------------|--|-----------------|-----------------------------|
| Software Free 30-day trial version | <table border="1" style="width: 100%; border-collapse: collapse;"> <tr> <td style="width: 50%;">How to DL and Install</td><td style="width: 50%;">How to DL and Install_T_E.pdf</td></tr> <tr> <td>Software(30Mb)</td><td>CellActivision_R1.02.01_64bit_demoInstaller.exe</td></tr> <tr> <td>Activation code</td><td>Req_ActivationCode_T_E.xlsx</td></tr> </table> | How to DL and Install | How to DL and Install_T_E.pdf | Software(30Mb) | CellActivision_R1.02.01_64bit_demoInstaller.exe | Activation code | Req_ActivationCode_T_E.xlsx |
| How to DL and Install | How to DL and Install_T_E.pdf | | | | | | |
| Software(30Mb) | CellActivision_R1.02.01_64bit_demoInstaller.exe | | | | | | |
| Activation code | Req_ActivationCode_T_E.xlsx | | | | | | |
| Software Licensed version | <table border="1" style="width: 100%; border-collapse: collapse;"> <tr> <td style="width: 50%;">How to DL and Install</td><td style="width: 50%;">How to DL and Install_L_E.pdf</td></tr> <tr> <td>Software(30Mb)</td><td>CellActivision_R1.02.01_64bit_unlimitedInstaller.exe</td></tr> <tr> <td>Activation code</td><td>Req_ActivationCode_L_E.xlsx</td></tr> </table> | How to DL and Install | How to DL and Install_L_E.pdf | Software(30Mb) | CellActivision_R1.02.01_64bit_unlimitedInstaller.exe | Activation code | Req_ActivationCode_L_E.xlsx |
| How to DL and Install | How to DL and Install_L_E.pdf | | | | | | |
| Software(30Mb) | CellActivision_R1.02.01_64bit_unlimitedInstaller.exe | | | | | | |
| Activation code | Req_ActivationCode_L_E.xlsx | | | | | | |

License version

- ② Start Installer in the folder, and install following the instruction wizard.




- ③ Start CellActivision after installation.



- ④ Window below appears.



- ⑤ Open the activation code request file,
"Req_ActivationCode_T_E.xlsx, which was downloaded at ①.

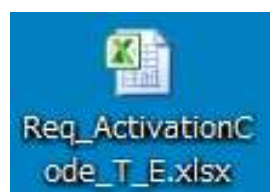

- ⑥ Transcribe “Site code” and “MID”.

CellActivision R1.01.01 Unlimited 64bit

Site code: DF4F26FD MID: D50E-1CCE-CF10-D3E9

CellActivision. Free 30-day trial version
Request for Activation code

Date: 2015/11/1

To: Yokogawa

Fill in the blank spaces and send this form to Yokogawa

| | | |
|-----------------|---------------------|--|
| Trial user name | Company | |
| | End-user name | |
| | e-mail | |
| Version | R1.01.01 | |
| Site code | MID | |
| DF4F26FD | D50E-1CCE-CF10-D3E9 | |

- ⑦ Transcribe the version number described in the file name of installer.

FH1-M3857LW.zip

File Edit View Tools Help

Organize Extract all files

File Name Type Compressed size

CellActivision_R1.01.01_64bit_unlimitedInstaller.exe Application 30,365 KB

CellActivision. Free 30-day trial version
Request for Activation code

Date: 2015/11/1

To: Yokogawa

Fill in the blank spaces and send this form to Yokogawa

| | | |
|-----------------|---------------------|--|
| Trial user name | Company | |
| | End-user name | |
| | e-mail | |
| Version | R1.01.01 | |
| Site code | MID | |
| DF4F26FD | D50E-1CCE-CF10-D3E9 | |

- ⑧ Fill in and save “Customer Information” column, and send the excel file “Req_ActivationCode_T_E.xlsx (Activation code request form)” to Yokogawa; “LSBioCAsupport@yg.jp.yokogawa.com”.

| CellActivision. Free 30-day trial version | | | |
|---|--|---------------------|--|
| Request for Activation code | | | |
| | | Date | |
| | | 2015/11/1 | |
| To: Yokogawa | | | |
| Fill in the blank spaces and send this form to Yokogawa | | | |
| Trial user name | | Company | |
| | | End-user name | |
| | | e-mail | |
| Version | | R0.01.01 | |
| Site code | | MED | |
| DF4F26FD | | D80E-100E-CF10-03E9 | |

- ⑨ Unlock CellActivision after receipt of “Activation code report” from Yokogawa; “LSBioCAsupport@yg.jp.yokogawa.com”, following the procedures explained below.

- ⑩ Start CellActivision.



- ⑪ Transcribe the “Activation code” described in the “Activation code report”, and click “Continue”.
CellActivision will be unlocked, and start.

| A | B | C | D |
|----|--|-------------------------------------|-----------|
| 1 | CellActivision. Free 30-day trial version | | |
| 2 | Activation code | | |
| 3 | | | Date |
| 4 | | | 2015/11/1 |
| 5 | To: Trial user | | |
| 6 | This is important information, please keep this form safe. | | |
| 7 | | | |
| 8 | | | |
| 9 | | | |
| 10 | | | |
| 11 | Trial user name | Company | 0 |
| 12 | | End-user name | 0 |
| 13 | | e-mail | 0 |
| 14 | | | |
| 15 | Version | R1.01.01 | |
| 16 | | | |
| 17 | Site code | MD | |
| 18 | PC version | D60E-10DE-0F10-03E9 | |
| 19 | Activation code | 020BF00B-FB3A3BDE-ACE7D64F-F40ECF79 | |
| 20 | | | |

CellActivision R1.01.01 Unlimited 64bit

Site code: 0F4F24F0 MD: D60E-10DE-0F10-03E9

Activation code:

☒ Unlock application

☐ Enter application

LOCKED

Cancel Continue >> English